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en Test kit for performing colorimetric tests on phosphate

Instructions for use: refer to the inside

This method can also be used for analyzing sea water after dilution (1+9).

Interferences: Large amounts of oxidizing reagents inhibit formation of the blue color complex. The following ions will not interfere: < 1 mg/L sulfide; < 10 mg/L citrate, heavy metals; < 50 mg/L Si.

Hazard warning/Disposal: Information can be found in the safety data sheet. You can download the safety data sheet from www.mn-net.com/SDS.

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DANGER Causes serious eye damage. | Wear protective gloves/eye protection. IF IN EYES: Rinse cautiously with water for several minutes. Remove contact lenses, if present and easy to do. Continue rinsing. Immediately call a POISON CENTER/doctor.

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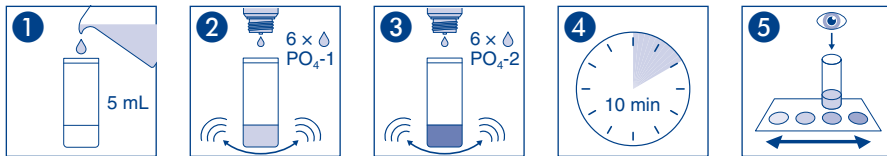
en Instructions for use

1. Rinse the test vessel several times with the water sample and fill to the ring mark (5 mL).
2. Add 6 drops of $\text{PO}_4\text{-1}$ and mix by swirling.
3. Add 6 drops of $\text{PO}_4\text{-2}$ and mix by swirling.
4. Wait 10 min.
5. Place the measuring vessel on the color chart and assign the value by comparison of the colour. Mid-values can be estimated.

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