
# IKA

designed for scientists

## A10 basic

Betriebsanleitung Ursprungssprache	DE	3			
Operating instructions	EN	14			
Mode d'emploi	FR	25			
Руководство по эксплуатации	RU	36			
使用说明	ZH	47			
取扱説明書	JA	58			
사용 설명서	KO	69			
คู่มือการใช้งาน	TH	80			
Indicaciones de seguridad	ES	91	Biztonsági utasítások	HU	101
Veiligheidsaanwijzingen	NL	92	Varnostna navodila	SL	102
Avvertenze per la sicurezza	IT	93	Bezpečnostné pokyny	SK	103
Säkerhetsanvisningar	SV	94	Ohutusjuhised	ET	104
Sikkerhedshenvisninger	DA	95	Drošības norādījumi	LV	105
Sikkerhetsinformasjon	NO	96	Saugos nurodymai	LT	106
Turvallisuusohjeet	FI	97	Инструкции за безопасност	BG	107
Instruções de segurança	PT	98	Indicații de siguranță	RO	108
Wskazówki bezpieczeństwa	PL	99	Υποδείξεις ασφαλείας	EL	109
Bezpečnostní upozornění	CS	100			

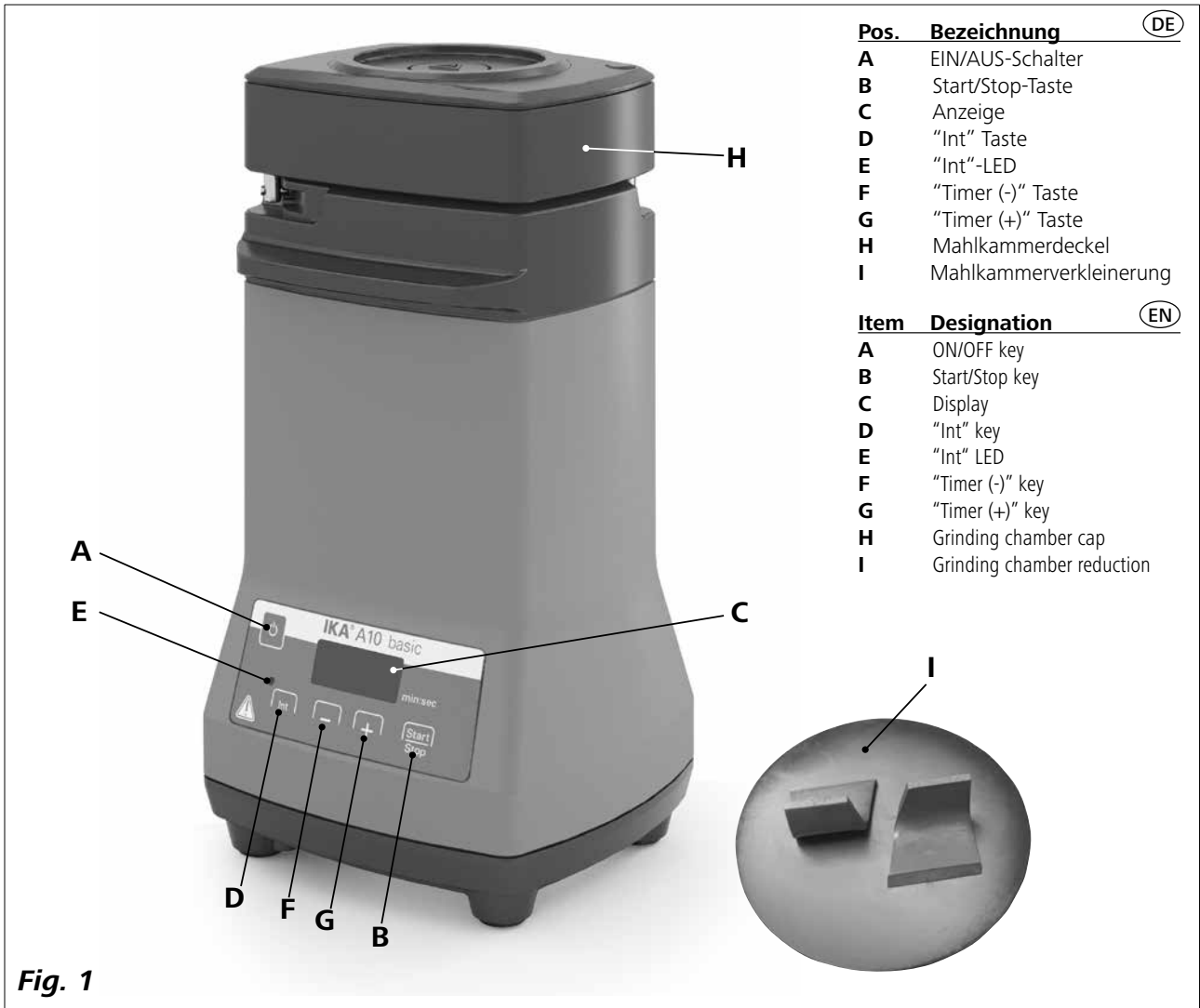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## Contents

	Page
Declaration of Conformity	14
Explication of warning symbols	14
Safety instructions	15
Correct use	16
Unpacking	16
Useful information	17
Commissioning	19
Timer function	19
Interval function	19
Exchanging and replacing the grinding tools	20
Maintenance and cleaning	21
Error codes	23
Accessories	23
Warranty	23
Technical data	24

## Declaration of Conformity

We declare under our sole responsibility that this product corresponds to the regulations 2014/35/EU, 2006/42/EC, 2014/30/EU and 2011/65/EU and conforms with the standards or normative: EN 61010-1, EN 6010-2-051, EN 61326-1, EN 60529 and EN ISO 12100.

A copy of the complete declaration of conformity or further declarations of conformity can be requested at [sales@ika.com](mailto:sales@ika.com).

## Explication of warning symbols

General hazard

This symbol identifies information **that is of vital importance for protecting your health and safety**. Disregarding this information may lead to health impairment and injuries.

This symbol identifies information **that is of importance for the technically correct functioning of the system**. Disregarding this information may result in damage to the instrument or to system components.

This symbol indicates information **which is important for proper use and ensuring that the operations of the instrument are performed efficiently and for using the instrument**. Failure to observe this information may result in inaccurate results.

## Safety Instructions

- **Read the operating instructions fully before starting up and follow the safety instructions.**
- Keep the operating instructions in a place where they can be accessed by everyone.
- Ensure that only trained staff work with the instrument.
- Follow the safety instructions, guidelines, occupational health and safety and accident prevention regulations.
- The instrument must be always supervised when in operation.
- Hearing protection must be worn when working with the instrument at high speed.
- Set up the instrument in a spacious area on an even, stable, clean, non-slip, dry and fireproof surface.
- The feet of the instrument must be clean and undamaged.
- The instrument is not suitable for hand held operation.
- The instrument may heat up when in use.
- Check the instrument and accessories for damage before each use. Do not use damaged components.

### DANGER

Please pay attention to avoiding hand injury when operate the instrument.

### DANGER

Always open the chamber carefully after use, as the material in the chamber may heat up due to transfer of energy during operation. Wear gloves if necessary or wait for material to cool down.

- Beware of the risk of flammable materials.
- **Do not** operate the instrument in explosive atmospheres, with hazardous substances or under water.
- The instrument doesn't start up again automatically following a cut in the power supply.
- Safe operation is only guaranteed with the accessories described in the "**Accessories**" chapter.
- Socket must be earthed (protective ground contact).
- The instrument could only be disconnected from the mains supply by pulling out the mains plug or the connector plug.
- The socket for the mains cord must be easily accessible.
- The voltage stated on the nameplate must correspond to the mains voltage.
- Protect the instrument and accessories from bumps and impacts.

### WARNING

Only process dry solid sample with the instrument. Never use it for processing liquid or wet samples.

- Do not crush explosive, toxic or noxious substances.
- Please observe the relevant safety instructions and guidelines regarding dust explosions.
- Please follow the relevant safety instructions and guidelines, and occupational health and safety regulations for use in the laboratory. When working with the mill, the user must wear his personal protective equipment.
- Always wear suitable protective equipment when embrittling the mill feed with coolants (dry ice).

### DANGER

Never cool the samples with liquid nitrogen inside **A10 basic** grinding chamber. If the samples have to be cooled with liquid nitrogen, please cool them first before filling into **A10 basic** grinding chamber.

- Ensure that the chamber cap is tightly closed.

### DANGER

Before removing the chamber cap, switch off the mill and wait until the motor stops. Beware of risk posed by trailing grinding tool.

- Always switch off the instrument and disconnect the plug before cleaning or replacing parts.

### DANGER

The grinding pot, grinding tool and rotor shaft may be hot after long grinding cycles. Allow the mill to cool down before cleaning or before changing grinding tool.

### DANGER

Beware of the sharp edged cutter in particular when changing grinding tool and when cleaning. Always wear your personal protective equipment (protective gloves, etc.).

- Ensure that pressure does not build up in the grinding chamber (particularly when using coolants).
- The instrument can only be opened by trained specialists, even during repairs. The machine is to be unplugged from the mains before opening. Live parts inside the machine may still be live for some time after unplugging from the mains.

## Correct use

### • Use

The **A10 basic** is a batch mill that allows for both impact milling and cutting milling.

#### Impact milling:

For use with hard, brittle material, dried material or material that is brittle because of cold (for example grain, coffee, etc.).

Here the grinding material is reduced by means of a beating rotor. The grinding material is thus broken up. The granularity of the final product is determined by the duration of the grinding, the level to which the container is filled and the physical properties of the material being processed.

#### Cutting milling:

For use with soft, fibrous grinding material (for example hay, paper, etc.).

Here the grinding material is cut by means of a rotating blade. In this case as well, the granularity of the final product is determined by the duration of the grinding, the level to which the container is filled and the physical properties of the material being processed.

**Intend use:** Tabletop device

### WARNING

The mill is not suitable for grinding wet material with too much water. Otherwise, the water may leak from the chamber and damage the device.

### • Range of use (indoor use only)

- Laboratories
- Pharmacies
- Schools
- Universities

This equipment is suitable for use in all areas except:

- Residential areas
- Areas that are connected directly to a low-voltage supply network that also supplies residential areas.

The safety of the user cannot be guaranteed:

- If the instrument is operated with accessories that are not supplied or recommended by the manufacturer
- If the instrument is operated improperly or contrary to the manufacturer's specifications
- If the instrument or the printed circuit board are modified by the third parties.

## Unpacking

### • Unpacking

- Please unpack the device carefully
- In the case of any damage a detailed report must be sent immediately (post, rail or forwarder).

### • Delivery scope

see **Fig. 2**

- **IKA A10 basic** (include **A10.1**) ①
- One **A10.4** grinding chamber reduction ②
- Two hose clips ③
- Two felt rings ④
- Tools kits ⑤
- Operating instructions ⑥
- A warranty card ⑦.

**Fig. 2**

## Useful information

The mill can be used in a wide range of applications. The list of grinding materials given below is not complete. It is possible to expand the range of applications further by means of grinding experiments with corresponding grinding procedures and sample preparations.

### **Impact milling:**

The **A10 basic** with a beater reduces soft, medium hard and brittle materials with a Mohs' scale of hardness of up to 5. For greater levels of hardness, the wear and tear on the pulverization tool is very high and grinding therefore becomes uneconomical. If the **A10 basic** is used for grinding materials with hardness greater than 5 Mohs, the **A10.3** hard metal beater, which is more resistant to wear and tear, must be used. This prevents the sample from being contaminated by bits of the grinding tool broken off by wear and tear.

Everything that breaks, is dry, and does not have a high fat content can be ground. Following is a list of substances that can be ground dry:

Wheat, barley, corn, malt, pectin, roasted coffee, nutshells, bones, ergot, peat, pharmaceutical, fertilizers, feed materials, spices, resin, potash, seeds, salts, cinders, tablets.

Tough grinding material must be cooled, for example by adding pulverized dry ice to the grinding pot.

The grinding material must be initially broken down into pieces with a feed size of 6 mm (the size of a grain of corn is ideal).

### **Cutting milling:**

The **A10 basic** with the **A10.2** star knife reduces bulky, elastic fibrous and soft materials with high cellulose content. Mixed goods such as trash must be free of iron and non-iron metals. The feed material must not be too moist or fatty. Otherwise the material may adhere to the sides of the grinding container.


Following is a list of some materials that can be reduced: Leaves, fibers, hops, cardboard, paper, hay, plastics, tobacco and roots.

The feed granularity should be no bigger than 10 mm.

**Note:** Materials that are not listed above should only be reduced in the **A10 basic** after confirmation from **IKA**, especially if there is uncertainty about the danger of explosion (dust explosion as a result of electrostatic charge).

### **Working with the A10.4 grinding chamber reduction:**

When the sample is very light or the sample quantity is very small (1- 5 ml), the user can insert the **A10.4** grinding chamber reduction into the grinding chamber for grinding as shown in **Fig. 3**.

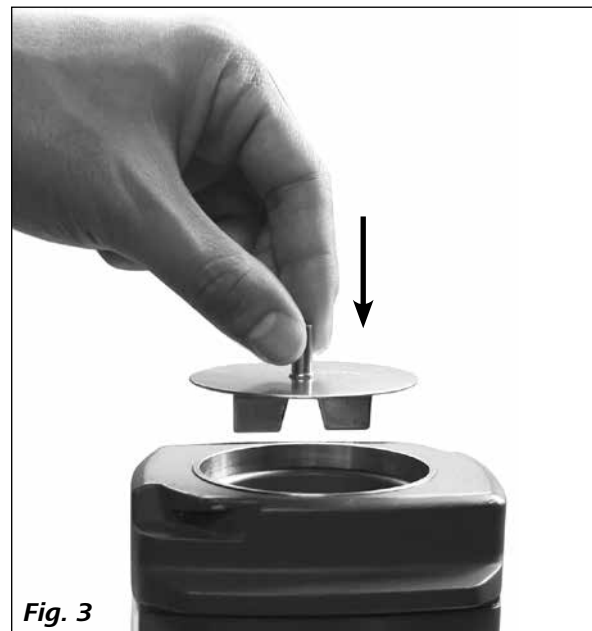**Fig. 3**

### **Working with cooling water:**

There are two hose couplings located at the bottom of the handle in order to connect the cooling water. Connect hoses as following:

- Remove the low part of the handle
- Connect the hoses to the couplings (we recommend hose  $\varnothing$  8 x 1,5 mm).
- Fix the hoses to the couplings with included clips.


**Fig. 4**

Then, tap water could flow into the device as cooling water from either hose.

**Note:** only water (max. 40 °C) could be used for coolant and water pressure should be less than 1 bar for the instrument. After the grinding, please ensure the cooling water is completely empty from the cooling chamber.

Cooling with tap water is usually sufficient for substances which become hot as a result of the grinding operation and thus are given to smearing or those that produce oil. The grinding chamber may only be filled halfway especially when samples containing oils and similar materials are ground.

**Note:** Moisture in material that is ground may evaporate because of the material temperature increasing. Then, the evaporated moisture will condense and flow out from the drain hole at the bottom of the instrument. This could prevent the motor from damage of moisture.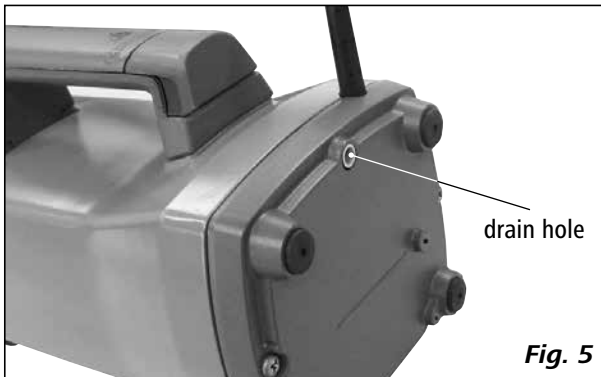

If more refrigeration is required, the ground matter should be cooled prior to being put into the grinding chamber.

Grinding dust may also leak out from this drain hole during grinding process.

### **Working with coolants – Dry ice cooling (CO<sub>2</sub>):**

Please follow the safety instructions.

Crushing causes the mill feed to heat up. This may not always be desirable as the mill feed can change when heated (e.g. increased oxidation, loss of moisture, evaporation of volatile components, etc.).

It may only be possible to crush mill feed which is viscous or has a high fat content by embrittling using coolants.

*The A10 basic offers following cooling processes:*

First put the feed product into the grinding pot and then add crushed dry ice.

Thoroughly mix the dry ice with the mill feed.


Only attach the chamber cap when the dry ice is completely vaporized, otherwise high pressure may build up in the grinding chamber.

Please note that the mill feed heats up again extremely quickly during grinding (it may reach room temperature again after approximate 10 seconds of grinding). You may have to repeat the process several times if the product is not adequately ground after one cooling process.

## Commissioning

Pay attention to the ambient conditions listed in the **"Technical data"**.

Check whether the voltage specified on the type plate matches the mains voltage available.

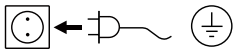The power socket used must be earthed (protective earth).

If these conditions are met, the device is ready to operate after plugging in the mains plug.

If these procedures are not followed, safe operation cannot be guaranteed and/or the equipment may be damaged.

### WARNING

Attach the grinding chamber cap by twisting it clockwise after filling grinding material. The cap is correctly attached only after **two "click"**!

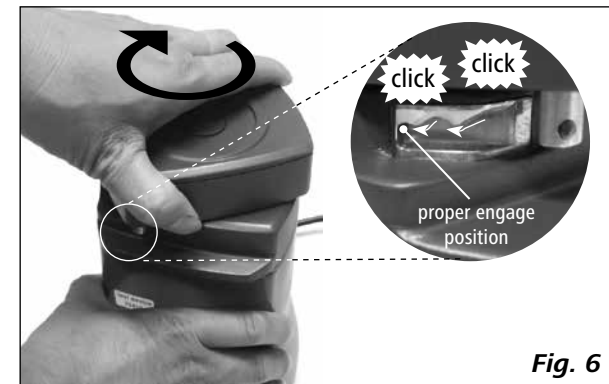
Fig. 6

When the mains plug is plugged in, all of the LEDs light up briefly during the self check.

Then, the software version is displayed.

After the software version, the right bottom LED lights up to indicate the power is connected.

Press ON/OFF switch (A), the display will show the **"Timer"** and **"Int"** setting.

Press the Start/Stop key (B), the instrument will start to run at the fixed speed 25000 rpm.

If the grinding chamber cap is not attached properly, the LED display three dashes (---) and grinding can not start.

Then, please check if the chamber cap is twisted to the stop position after two **"click"**.

If three dashes (---) don't disappear after the cap is correctly attached, please send the instrument for repair.

**Note:** When the interval function is activated by pressing **"Int"** button (D), the **"Int"** LED (E) lights up.

## Timer function

In stop status, press the Timer **"-"** button (F) and **"+"** button (G) to adjust the timer value. Adjusting magnitude of the timer is 1 second.

Keep pressing the Timer **"-"** button (F) or **"+"** button (G), the time will change continuously. The maximum timer value is 5 minutes (5:00).

When the timer is activated, the value on the display will countdown.

When the timer is deactivated, the instrument will automatically count from 0 second (0:00) to 5 minutes (5:00) and then the unit will stop running.

### WARNING

After the instrument runs 5 minutes, it should stop 10 minutes for cooling the motor. Then the user could start the instrument again by pressing Start/Stop button (B) again as request.


## Interval function

In stop status, press the button **"Int"** key (D), the interval function will be activated. The **"Int"** LED (E) lights up.

If timer function is activated, when the interval function is activated, the instrument run 15 seconds and stop 1 second repeatedly until the setting time run out.


If timer function is deactivated, when the interval function is activated, the instrument run 15 seconds and stop 1 second repeatedly until 5 minutes (5:00) elapse.

## Exchanging and replacing the grinding tools

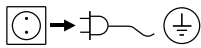 Disconnect the mains plug prior to exchanging or replacing the grinding tools!


Exchanging and replacing the grinding tools as following: Hold the grinding tool securely in place with the bracket. Loosen the hexagonal retaining nut of the grinding tool with socket wrench **SW 8** as indicated in following **Fig. 7**.

**Fig. 7**
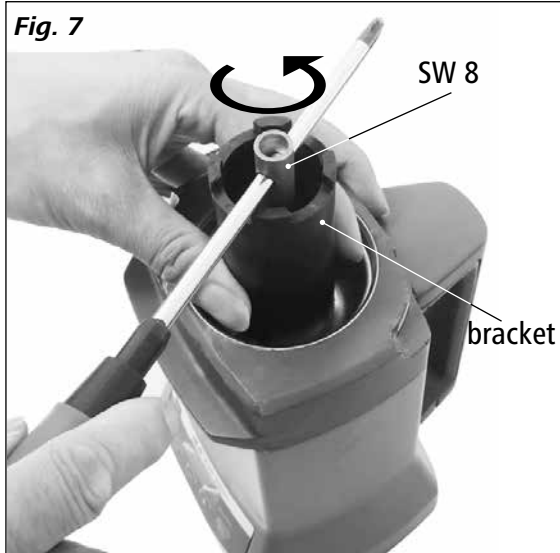


Then, the hexagonal nut and the grinding tool can be removed from the chamber.

**Fig. 8**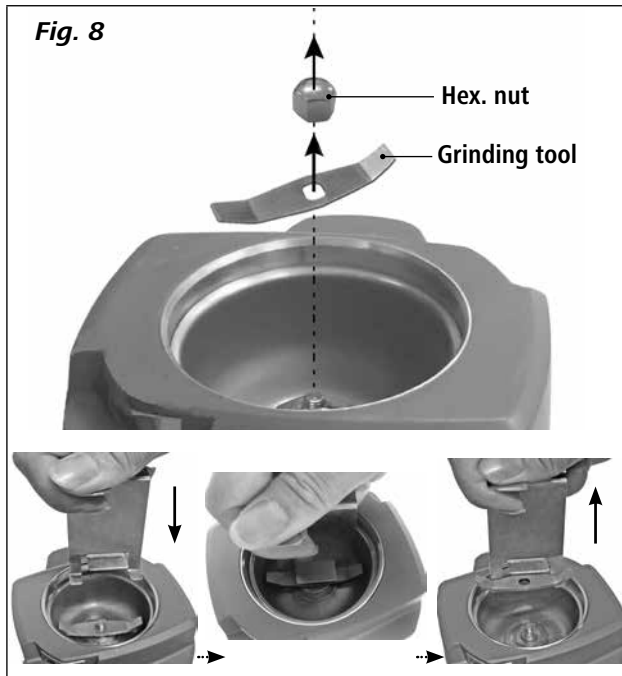

**Note:** When removing the **A10.2** Star-shaped beater, use the hooks on the bracket to hold the cutter. Then, the cutter can be taken out of the chamber (see **Fig. 9**, **Fig. 10**).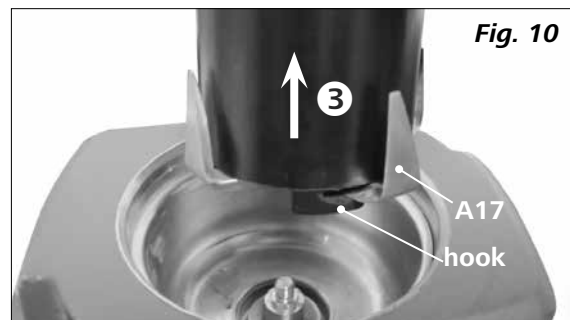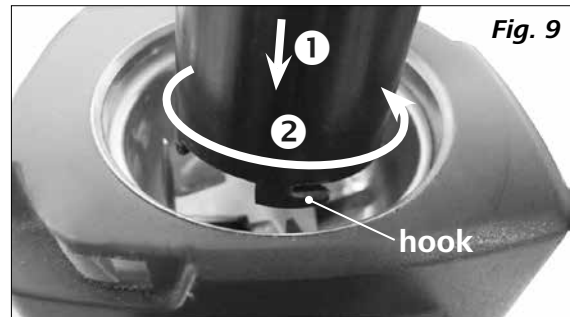

 **DANGER**

**The edges of the blade are sharp. It should only be handled with protective gloves. Please observe the safety instructions.**

After set the new or another grinding tool in place, tighten the hexagonal nut in clock wise direction. To do this, use the tools as shown in **Fig. 7**.

 **DANGER**

Note that when changing the tool directly after a grinding procedure, the pulverization tools are hot. The mill must first be allowed to cool off.


Before assembling a new or another pulverization tool, make certain that all parts are clean.

## Maintenance and cleaning

### Maintenance

If grinding tools wear out over the course of time, especially as a result of constant grinding of very hard materials, so that rapid grinding with the appropriate quality is no longer possible. The worn out parts must be replaced. For replacing grinding tools, see section “**Exchanging and replacing the grinding tools**”.

### Cleaning

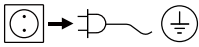
 Disconnect the mains plug prior to cleaning!

Dirty parts in the pulverization area can be cleaned with a cleaning brush or dry cloth.


If there is heavy accumulation of dirt, the grinding tool and grinding pot can be disassembled and cleaned in a dishwasher.

Disassemble the grinding tool and container as following:

- Disassemble the grinding tool (see “**Exchanging and replacing the grinding tools**” section).
- Disassemble the grinding pot with socket wrench **SW 17**.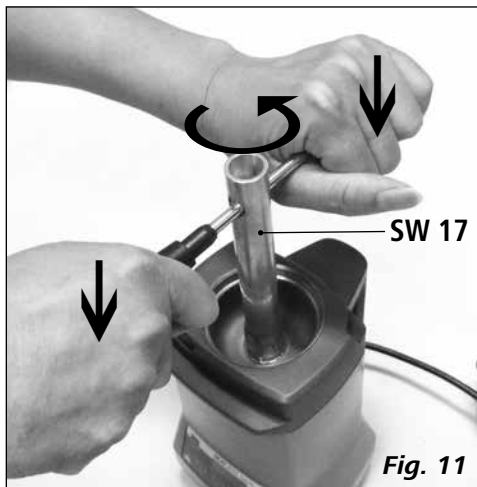

Then, grinding pot and grinding tool can be disassembled and cleaned with a dishwasher or by hand.

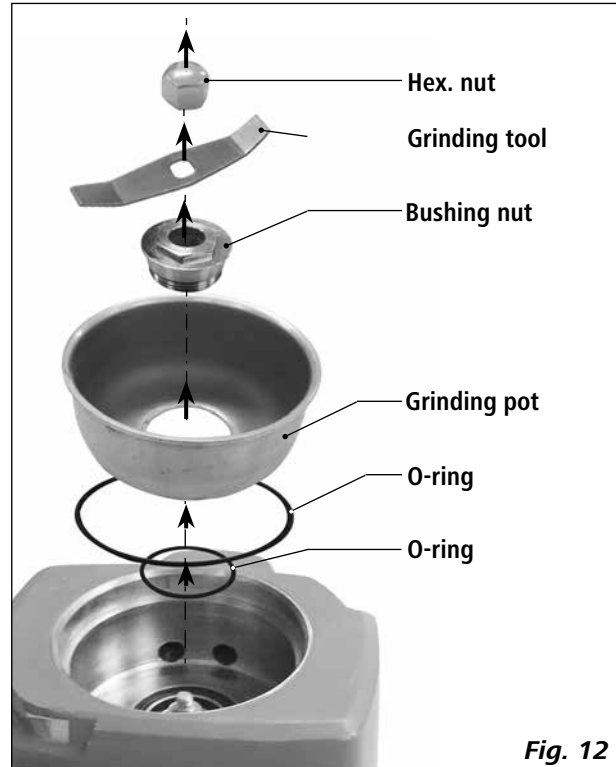**Note:** The chamber cap cannot be cleaned in a dishwasher.

After cleaning, all parts can be assembled in reverse sequence.

**Note:** Please pay attention to O-rings during all process. When O-rings wears out, replace them timely.

For dust-free and avoid cross contamination, we recommend to replace the felt ring in the bushing nut after every grinding with help of a small screw driver.

If grinding dust is found in the cavity below the grinding container, please clean it with included cleaning brush (see **Fig. 14**).

**Fig. 14**

You can also clean it with a vacuum cleaner (see **Fig. 15**).

**Fig. 15**

Interior of the device can be cleaned with compressed air via the drain hole (see **Fig. 16**).

**Note:** be aware of the threat by dusting!

**Fig. 16**

If there is heavy accumulation of dirt on this area, please wiped clean with a moist cloth. Never clean it in water or in a dishwasher.

Use only cleaning agents which have been approved by **IKA** to clean the instrument and all parts.

**Dirt**

- Dyes
- Construction materials
- Cosmetics
- Foodstuffs
- Fuels

**Cleaning agent**

- isopropyl alcohol
- water containing tenside/  
isopropyl alcohol
- water containing tenside/  
isopropyl alcohol
- water containing tenside
- water containing tenside

For materials which are not listed, please request information from **IKA**.

Wear protective gloves during cleaning the instrument. Electrical devices may not be placed in the cleansing agent for the purpose of cleaning.

Do not allow moisture to get into the equipment when cleaning.

Before using another than the recommended method for cleaning or decontamination, the user must ascertain with **IKA** that this method does not destroy the instrument.

**Spare parts order**

When ordering spare parts, please give:

- machine type
- serial number, see type plate
- item number and designation of the spare part, see [www.ika.com](http://www.ika.com), spare parts diagram and spare parts list,
- Software version.

**Repair**

**Please send equipment for repair only after it has been cleaned and is free from any materials which may constitute a health hazard.**

For this, you should request the “**Decontamination Certificate**” from **IKA**, or use the download printout of it from the **IKA** website [www.ika.com](http://www.ika.com).

If you require servicing, return the equipment in its original packaging. Storage packaging is not sufficient. Please also use suitable transport packaging.

## Error codes

The fault is shown by an error code in the LED display (C) as following if the error occurs.

Proceed as follows in such cases:

- ☞ Switch off the device using the main switch
- ☞ Carry out corrective measures
- ☞ Restart the instrument

Error code	Cause	Effect	Solution
---	The cap is not closed properly	instrument does not start	- Close the cover properly
<b>E3</b>	Temperature inside instrument is too high	motor off	- Switch off the instrument and allow it to cool down
<b>E4</b>	Motor blockage or overload	motor off	- Switch off the instrument - Decrease the motor's load and reduce the material to be ground
<b>E25</b>	Safety relay fault	motor off	- Send the instrument for repair
<b>E86</b>	Voltage is too low	motor off	- Switch off the instrument - Check the Mains voltage

If the actions described fails to resolve the fault or another error code is displayed then take one of the following steps:

- Contact the service department
- Send the instrument for repair, including a short description of the fault.

## Accessories

- A10.1** Standard beater
- A10.3** Hard metal cutter

- A10.2** Star-shaped knife
- A10.4** Grinding chamber reduction

## Warranty

In accordance with **IKA** warranty conditions, the warranty period is 24 months. For claims under the warranty please contact your local dealer. You may also send the machine direct to our factory, enclosing the delivery invoice and giving reasons for the claim. You will be liable for freight costs.

The warranty does not cover worn out parts, nor does it apply to faults resulting from improper use, insufficient care or maintenance not carried out in accordance with the instructions in this operating manual.

## Technical data

Operating voltage	<b>VAC</b>	220–240 ± 10 % 100–120 ± 10 %
Frequency	<b>Hz</b>	50 / 60
Motor rating input	<b>W</b>	300
Motor rating output	<b>W</b>	240
Rated speed	<b>rpm</b>	25000 ± 2 %
Maximum circumferential speed	<b>m/s</b>	73
Motor		brushless motor
Maximum useful volume	<b>ml</b>	50
Maximum feed hardness	<b>Mohs</b>	5
Maximum feed grain size	<b>mm</b>	6 – impact milling 10 – cutting milling
Parts in contact with product		stainless steel
Short time operation (KB)	<b>min</b>	5 / 10 (5 min. ON / 10 min. OFF)
Cooling water pressure	<b>bar</b>	max. 1
Cooling water temperature	<b>°C</b>	max. 40
Timer max.	<b>min</b>	5
Time setting accuracy	<b>sec</b>	1
Interval	<b>sec</b>	15 / 1 (15 seconds ON / 1 second OFF)
Protection class according to EN 60529		IP 41
Ambient temperature	<b>°C</b>	+ 5 ... + 40
Ambient humidity (relative)	<b>%</b>	80
Operation at a terrestrial altitude	<b>m</b>	max. 2000
Dimension (W x D x H)	<b>mm</b>	130 x 145 x 250
Weight	<b>kg</b>	2,9

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**IKA Works (Thailand) Co. Ltd.**
**IKA Works Kenya Ltd.**
**IKA Works Kampala Limited**